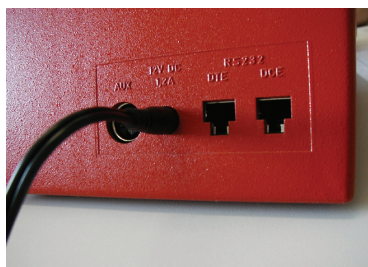
Guida rapida al FOODLAB . wineLab

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
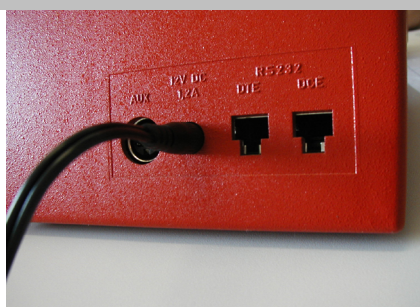
FOODLAB and WineLab short guide

To switch on the instrument insert the plug...

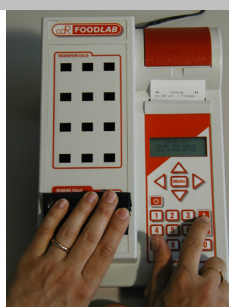... and press "ON" button.

DISPLAY
Calibrate Cover the wells and press ENTER

The instrument may request automatic calibration. Cover the reading cells with the pad provided and press ENTER. Wait for some seconds.

The result of the check is shown on the display

DISPLAY
Calibrate
2 ok 4 ok 6 ok
1 ok 3 ok 5 ok

On FOODLAB Fat the channel 2 and 5 are --
On WineLab the channel 2 and 6 are --
Press ENTER. The instrument is operative.
If the calibration fails, the system will indicate "KO" beside the well that has encountered the problem.
Please contact CDR s.r.l. for assistance.

DISPLAY
Analysis
test n°1
test n°2
...

Press the "0" key to move from the screen showing the wells' status (stand-by) to the analysis list. Press again "0" to go back to the stand-by screen.

After an analysis session, to start a new analysis and cancel the results press ENTER...

... and then the ▽ arrow.
Press "0" if you want to go back to the stand-by screen and select an analysis that works with a different well.

The ▽ arrow is used to cancel or abort an action or go back to previous step.

DISPLAY
Menu
Edit...
Setup...
Autostart

To access the settings MENU press in rapid sequence the following keys "ENTER" "1" "2" "3" "ENTER". Then press MENU.
In WineLab press in rapid sequence the following keys "0" "1" "2" "3" "ENTER". Then press MENU.

CORRECT PROCEDURE**INCORRECT PROCEDURE**

Insertion of a new tip: push the piston with force right to the end of its travel so that the tip adheres well.

Do not leave a space between the tip and the pipette: push it fully in.

Insert the tip cover with the fingers and push fully home until a click is heard.

Do not leave space between the tip cover and the pipette.

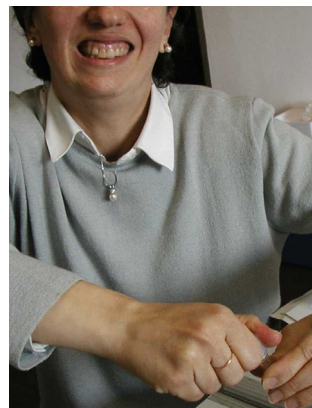
Turn the piston until the required volume is centred by the arrow (e.g. 5 μ l = 050).

Before making an analysis, check that the volume in the pipette is correct.

CORRECT PROCEDURE**INCORRECT PROCEDURE**

The caps must be opened carefully by putting pressure on the right and then left. Do not touch the internal parts (soaked) of the cap with the hands.

Do not open it by using force: it will never open.

When taking the sample, always agitate by inverting the container with the sample and immerse the tip keeping the piston lowered.

Do not raise the piston in jerks but take the sample SLOWLY.

The cleanliness of the tip is fundamental for making a correct analysis: clean it by sliding the tip on paper.

Do not touch the point of the tip so as not to take the sample from inside.