

Updating your CDR analyzer

CDR FoodLab®, CDR WineLab®, CDR BeerLab®, CDR CiderLab, CDR OxiTester, CDR KombuchaLab, CDR PalmOil Tester

For 222 Model with software version 3.0 or later.

For 242 Junior model with software version 3.6 or later.

During the updating procedure you will be invited to register your analyzer if not already done.

1. Turn analyzer on and allow it to warm up




2. Select "Settings"




3. Select "Update"



4. Follow prompt on screen and plug your analyzer to a computer using the provided white usb cord

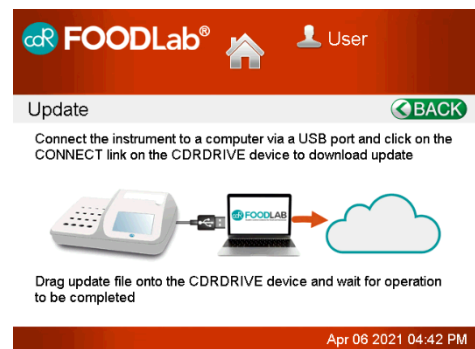


5. Open the CDRDRIVE folder on your computer

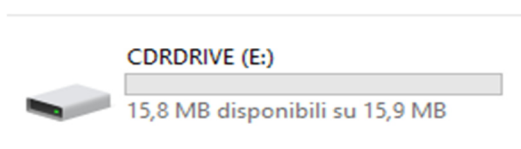


6. Click the .connect link in the CDR DRIVE

You will be routed to CDR's resource website

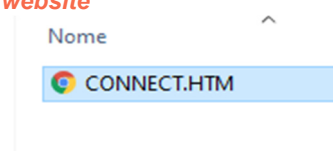

7. At the top of your screen, click on “Your instrument is not updated, please update it!”

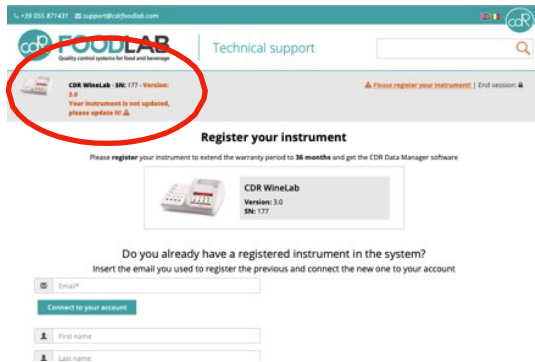

8. Click to download the latest update

Software updates




Software version

CHANGE LOG:


- new analysis available (opt)
- resolution of various bugs
- minor changes

Follow the steps below to update

9. Drag downloaded file into the CDRDRIVE folder in order to begin the download

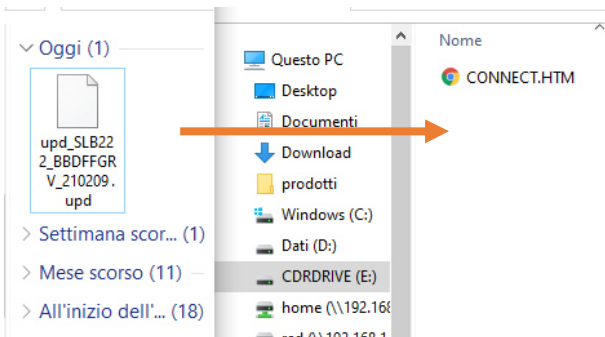

10. Wait the updating procedure is completed. At the end the analyzer reboots

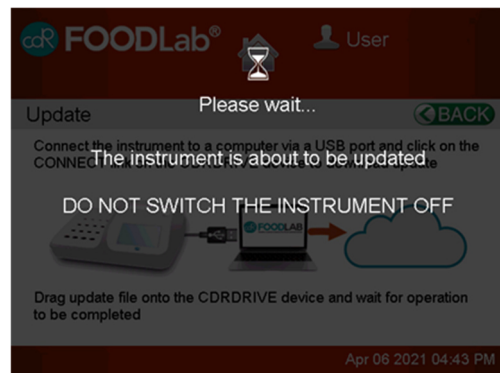

11. Disconnect the analyzer from PC and connect it again. In CDRDRIVE folder click on connect once again

